

NEW



SITVue™

DAB Detection System

SITVue™/DAB Detection System is a powerful Intensification system that significantly enhances chromogenic signals. This system can be integrated into standard immunohistochemistry (IHC) staining methods.


SITVue™/DAB Detection System is an enzyme-mediated reaction that utilizes horseradish peroxidase (HRP) to catalyze the deposition of two separate Linkers applied sequentially onto tissue sections or cell preparation. The deposited Linkers can be detected with a streptavidin-peroxidase conjugate followed by a reaction with a peroxidase substrate/chromogen solution such as diaminobenzidine (DAB).

The SITVue™/DAB Detection System results in a significant increase in sensitivity compared to standard IHC detection methods, while maintaining similar specificity.




- FAST**
Results in 30min
- DEPENDABLE/CONSISTENCY**
Optimized for use with Diagnostic BioSystems Concentrated or Ready to Use IHC antibodies
- RELIABLE**
Significant increase in staining intensity

SITVue™ /DAB Detection System

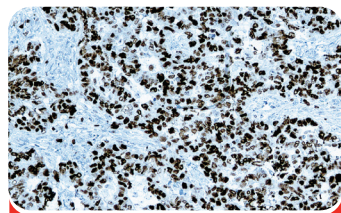


Ki67

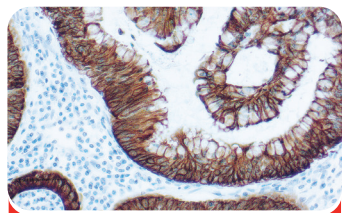


Cytokeratin Pan

Two Steps Detection System

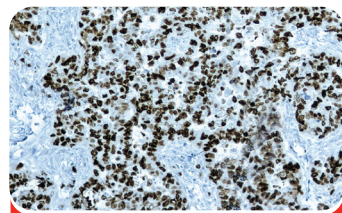

Ki67

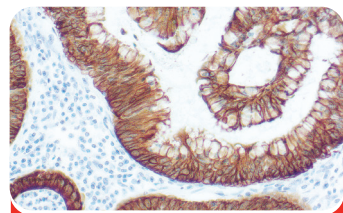

Cytokeratin Pan

Ordering Information

#	Cat. No	Volume
1	SIT-1000D	1000 Tests
2	SIT-100D	100 Tests

Step	Staining Procedure	Two Steps Detection system Incubation Time	SITVue™ /DAB Detection System Incubation Time
Peroxidase Block	A. Incubate slides in Peroxidase Block.	5 min.	1min.
	B. Rinse slides with Immuno Wash Buffer three (3) times, for 1 min. each time.	3 x 1 min.	
Pre-Blocking (optional)	A. Add 2 drops (100 µL) or enough volume of Pre-Blocking Solution to cover the tissue section.	10 min.	N/A
	B. Drain or blot off solution. Do not rinse.		
Primary Antibody	A. Incubate with Primary Antibody, prepared according to the manufacturer's recommended protocol at the desired concentration. Concentrated Primary Antibodies may be diluted using Primary Antibody Diluent.	30 – 60 min.	5 min.
	B. Wash slides with 3 changes of Immuno Wash Buffer.	3 x 1 min.	3 x 10 second.
Linker 1	A. Apply the Linker 1 and incubate.	10 min.	5 min.
	B. Rinse slides with 3 changes of Immuno Wash Buffer.	3 x 1 min.	3 x 10 second.
Linker 2	A. Incubate the tissue with Linker 2.	N/A	5 min.
	B. Wash slides with 3 changes of Immuno Wash Buffer.		3 x 10 second.
Tracer	A. Incubate the tissue with Tracer.	10 min.	5 min.
	B. Wash slides with 3 changes of Immuno Wash Buffer.	3 x 1 min.	3 x 10 second.
Stable DAB/ Plus	A. Prepare the Stable DAB/Plus substrate working solution (see above).		
	B. Incubate tissue with prepared Stable DAB/Plus substrate solution. Monitor level of staining to determine optimal time of incubation.	5 – 10 min.	5 min.
	C. Rinse slides with 3 changes of water.	3 x 1 min.	3 x 10 second.
Counterstain	A. Incubate tissue with Counterstain (Hematoxylin K097), according to manufacturer's recommendation or standard laboratory protocol.	~1 min.	~1 min.
	B. Wash slides with water 3 times, followed by 1 time in Immuno Wash Buffer, then 1 time in water.	3 x 1 min. H2O 1 x 1 min. Buffer 1 x 1 min. H2O	3 x 10 second. H2O 1 x 10 second. Buffer 1 x 10second. H2O
Total staining time		90-125 min	< 30 min
Dehydrate & Coverslip	A. Dehydrate tissues through graded ethanol series, followed by xylene series. B. Apply coverslips with permanent mounting medium.		



United States (Headquarters)

6616 Owens Drive, Pleasanton, CA 94588, USA
Toll Free: (888) 896-3350, customersupport@dbiosys.com



Contact your distributor for more information and visit us at www.dbiosys.com

Europe

Germany
Lake Constance
Cell: +49 7557 929 3915
wolfgang.vogel@dbiosys.com

Mexico

Avenida Aztecas # 95, Casa 4
Colonia Pueblo de los Reyes
Alcaldia Coyoacan
CDMX, C.P. 04330
Cell: +525535709693
a.herrera@dbiosys.com

India

612, Eden Square
St. John's Road
Secunderabad – 500003
Cell: +91 9958293222
anandam@dbiosys.com

Canada

32 Lemsford Drive
Markham, ON L3S 4H4
Cell: +1 416 219 2035
hosna.mujadidi@dbiosys.com